

<b>Title:</b>	<b>BCR Commands</b>	<b>ID:</b>	
		0011	
<b>Date in:</b>	<b>Response:</b>	<b>Model:</b>	<b>Author:</b>
2002-06-22	2002-06-22	-	CMa

BCR STX,LPX

On this carrousel you position the carrousel and lift to the plate you want to read:

"WR DM0 x"  
"WR DM5 y"  
"ST 1910"

This will position the lift to the plate xy. When you're there

RD 1915 -> '1'

you may read the barcode by sending

"LON"+ cr (9600 Bd, even, 7, 1 on 2nd comport).

Once the 1910 is set you can go to other locations by simply adjusting the value in DM5 or DM0.

**Some other commands for the BCR:**

"LOFF" (laser off)  
"MOTOROFF"  
"MOTORON"  
"TEST1" (continuous reading)