

Title:	How to detect CO2 option	ID:	
		0132	
Date in:	Response:	Model:	Author:
2004-05-19	2004-05-19	-	CMa

Q:

We currently have a Liconic 42 pos. incubator part of one of our demo systems.

We would like to be sure that the CO2 regulation is available and already fully installed on this equipment.

Could you please tell me how could we check it ?

(We do not use CO2 for demo but this unit could be sold to a customer soon and we need to be sure of its functionalities).

A:

If there is a third display you have CO2. I would fill some CO2 through the hole in the front glass door (~10 seconds) and watch the CO2 reading to rise.