

Title:	Humidification Control	ID:	
		0141	
Date in:	Response:	Model:	Author:
2004-06-24	2004-06-24	-	CMa

Q:

Yet another question:

What does this RH option on the StoreX give us?

Is this an active humidification system or just a monitor?

With this option do you still use a water pan at the bottom of the incubator?

A:

Yes, it is active controlled but it will not allow a free choice of the humidity level. This means that it will always be very humid but it will increase humidification on a door or gate access in order to accelerate recovery. The water pan is still there.