

<b>Title:</b>	<b>UPS STR240</b>	<b>ID:</b>	
			0284
<b>Date in:</b>	<b>Response:</b>	<b>Model:</b>	<b>Author:</b>
2007-01-23	2007-01-23	STR240	CMa

## Q:

I need your help regarding one of our customers that is currently using the STR240 incubator integrated to one of our BioCel. They would like to connect the incubator to an UPS in order to have the incubator running if they face an electrical shortcut. Could you please let me know which UPS you recommend for the STR240?

## A:

We have no recommendation for a specific UPS. However, V11 splits high power (climate) and low power (handling). The power requirements for the handling is 50..100 Watt. For the high power you will need ~1000W peak. The power consumption depends a lot on the climate settings. To give you an idea I added some plots taken from an STX220.