

| | | | |
|-----------------|------------------------|---------------|----------------|
| Title: | Swap Commands 2 | ID: | |
| | | 0229 | |
| Date in: | Response: | Model: | Author: |
| 2003-07-15 | 2003-07-15 | - | CMa |

Q:

We've got an Overlord installation which uses a Liconic carousel with a 'Turn' Station. The driver we have written does not use the OCX, its an old driver which uses the serial commands. Do you have a command that will allow me to activate the Turn Station?

A:

I'm assuming that you mean a Swap Station. Then Turn Station (Landscape/Portrait) is activated automatically on ever access of the internal robot. The Swap Commands are:

ST 1912: Swap Out

RS 1912: Swap In

RD 1912: Read Swap Status

The Flags to check for a plate are

RD 1813 Read Sensor on Nest 1

RD 1807 Read Sensor on Nest 2

Note that the Swap Station operates in Slave Mode. This means that the surrounding systems decides what Nest is in or out.