

Title:	Heaters do not Work 1700 Flag	ID:	
		0316	
Date in:	Response:	Model:	Author:
2008-10-28	2008-10-28	-	CMa

Q:

Thanks to your advise to send "RS 1700", 1700 is now "0", PLC 513 is ON and switched heater gets AC100V. Temperature in the chamber is rising toward the set point.

Thank you very much for your support.

Looking back what happened, both PLC taken from Steri-CultR in stock and a new PLC originally had "1700=1" setting. Our engineers now gain knowledge to watch "1700" when PLC is exchanged.

Is there any other parameter (flag) they have to be careful when PLC is exchanged ?

A:

Yes, you have to be very careful when going into this mode. In this mode the outputs are no longer controlled by the internal logic. This means that if an output is set it remains set until you reset it (or you leave this mode). It is planned to remove this feature since I consider it to be too risky.