

Title:	Adjusting CO2	ID:	
		0324	
Date in:	Response:	Model:	Author:
2007-08-07	2007-08-07	-	RKi

Q:

The CO2 is showing a wrong measurement when compared to a reference measurement.

A:

First check the CO2 reading at a low level like 2%.

- 1) If the display shows the two percent and the reference CO2-Gauge shows a value slightly above the Display can be adjusted.
- 2) If the measurement of the reference Gauge shows a huge difference, like the real value in the chamber is at 10% or above, a component of the CO2 conditioner might be defect

In Case 1) Follow the Instructions show in the Q&A for Adjusting the CO2

In Case 2) Check Sensor if still working. Check filter if water is in it.