

Title:	Shovel Touches Plates in Cassette			ID:
				0338
Date in:	Response:	Model:	Author:	
2008-01-08	2008-01-08	-	CMa	

Q:

carousel Model no: 189MP163DWP
REF 10612312
S/N 611000001

When shovel are picking plates from the towers the shovel are very close to bottom of the plates in some positions.

Some time the shovel touch the bottom of the mtp which result in not picking up the plate correctly.

How do I adjust the space between the shovel and the bottom of mtp in the towers ?

A:

There is a set screw for the lift wagon at the base plate next to the lift column. If you turn this screw in, the lift will initialize lower. There is also the possibility that the shovel is bent in which case the shovel is not parallel to the MTP