


| | | | | |
|-----------------|------------------------------------|---------------|----------------|------------|
| Title: | RH Crosstalk on CO2 Readout | | | ID: |
| | | | | 0389 |
| Date in: | Response: | Model: | Author: | |
| 2011-10-27 | 2011-10-27 | | CMA | |

The issue between RH and CO2 going to 10+ percent. Unit has steam generator. Please check.

Problem not know. I would first measure the signal on the CO2 sensor (5.0 .. 6.0 VDC) to see if originates from the sensor.

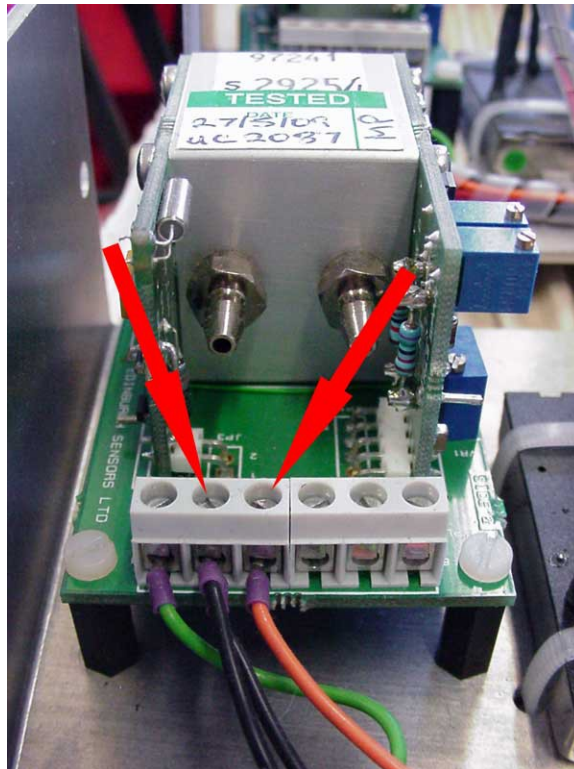

If not, take measurement at the display next.