

Title:	DF Defrost Procedure			ID:
				0077
Date in:	Response:	Model:	Author:	
2006-07-07	2006-07-07	STX1000	CMa	

Complete Defrost Procedure STX1000-DF Deep Freezer



Make sure to place container under drain output. "ST 1405" command starts complete defrost cycle. When defrost is complete this will also turn off the compressor.



The actual evaporator temperature is displayed on the 3rd display After ~5..10 min. water starts flowing. At the end of the defrost cycle the water flow should




slow down or should stop. The end of the defrost cycle is reached when the evaporator temperature is larger then 20°C.

To heat up unit remove one or two cassettes and place electrical heater on carrousel. The heater power should be 1..2kW. Ideally the heater has a built-in



Temperature regulator. Close door after placing heater. To avoid excessive frost inside the chamber do not leave door open. It is important to close door as quickly as possible.

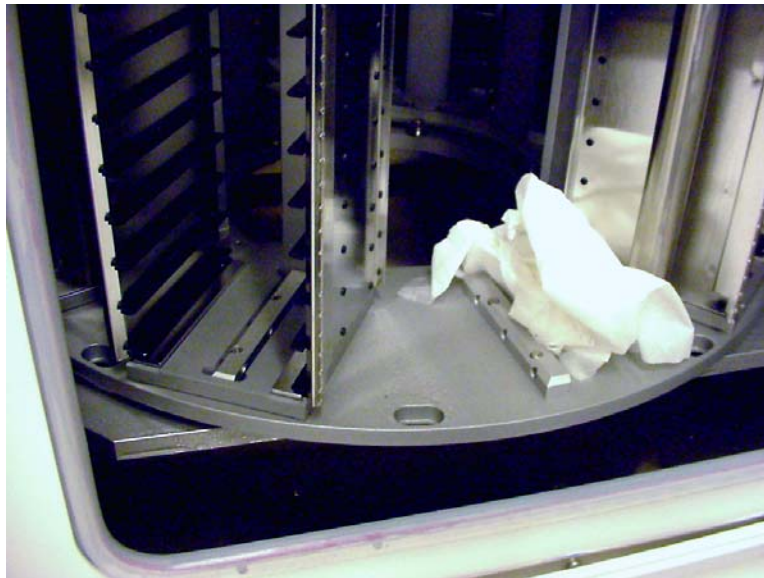

After ~15..30 min. interior reaches room temperature. Dry condense water using a soft cloth.