


Title:	Alarm Connector	ID:	
		0102	
Date in:	Response:	Model:	Author:
2010-09-30	2010-09-30	-	CMa

Alarm Connector User Information



Location of Alarm Socket on Connector Panel

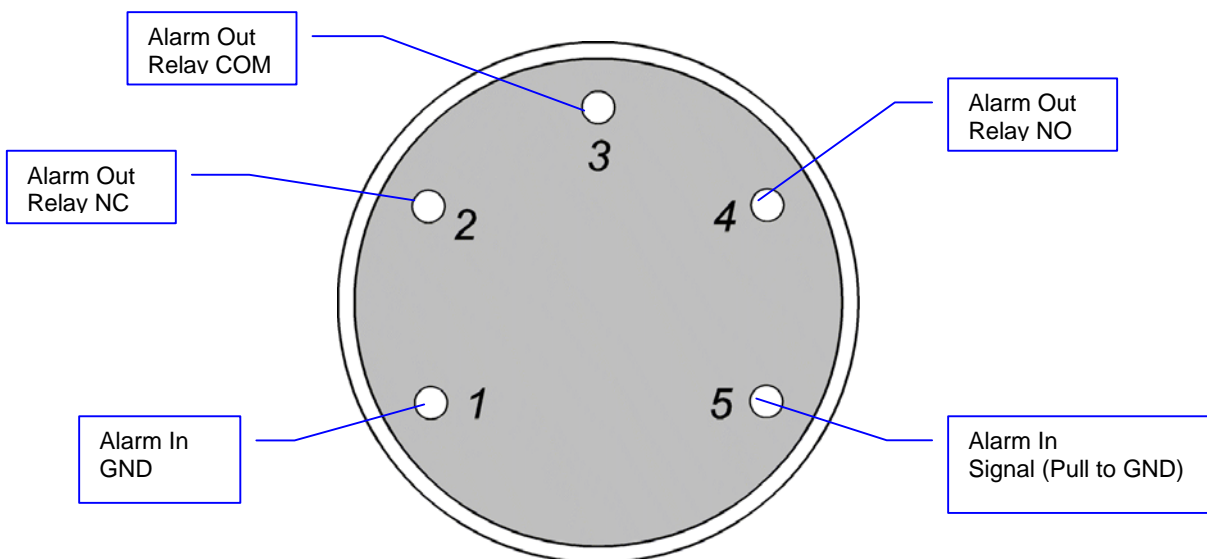

Alarm Socket Pin Layout seen from Connector Side

Connector Type Information

Connector Type Information: DIN Connector 5-Pin / 270°



Connector Side View



Connector Pin View

Low Level Alarm Paramter Settings

- Alarm Output for Temperature out of band is 600
- Alarm Set point DM900, value: $(t+500)*32$ [1/10°C]
- Band for temperature is DM130 [1/10 °C]
- Delay for Temp is T130 [1h]
- Alarm turns on after T130 time [1/10s]. As soon as temp. is within band.
- Reset Alarm: "RS T130" resets alarm for T130 (1 hour)