


Title: Current Resistors Stepper Card			ID:
			137
26.08.14	Response:	Model:	Author:
		-	CMA

Current Resistors Stepper Card

Stepper Motor Controller Cards NMC V1.xx

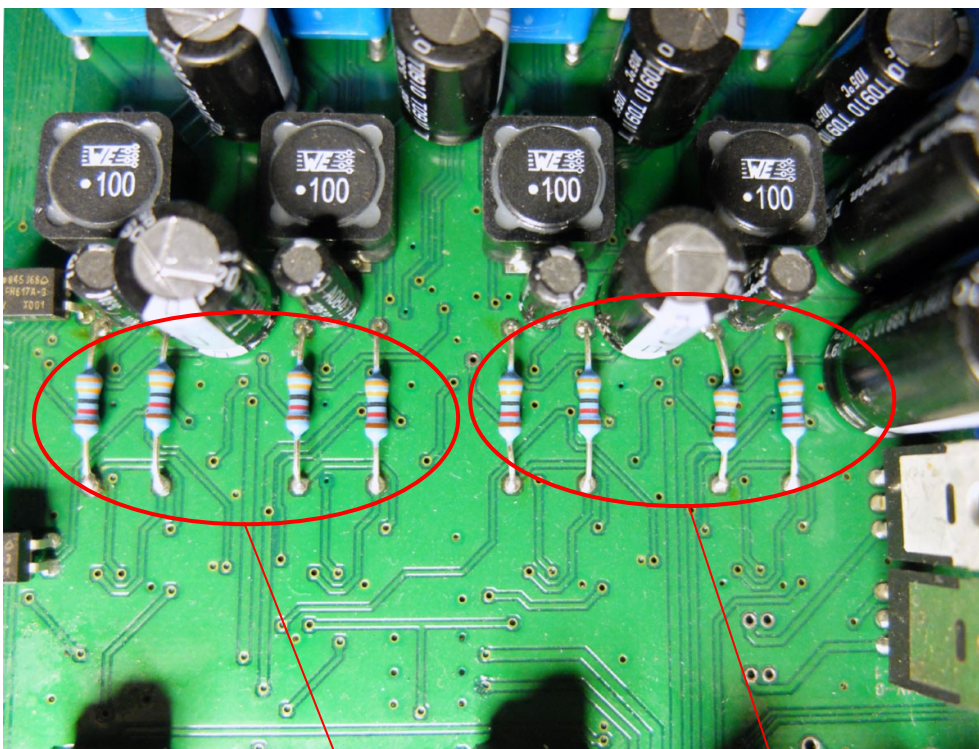


Current Resistors
(4) Channel 2

Current Resistors
(4) Channel 1

Resistor	Current
8.2k	1.1A
10k	1.4A
15k	2.1A
22k	3.2A
33k	4.5A

Resistors must be matched / paired within each axis!

Stepper Motor Controller Cards NMC V2.xx

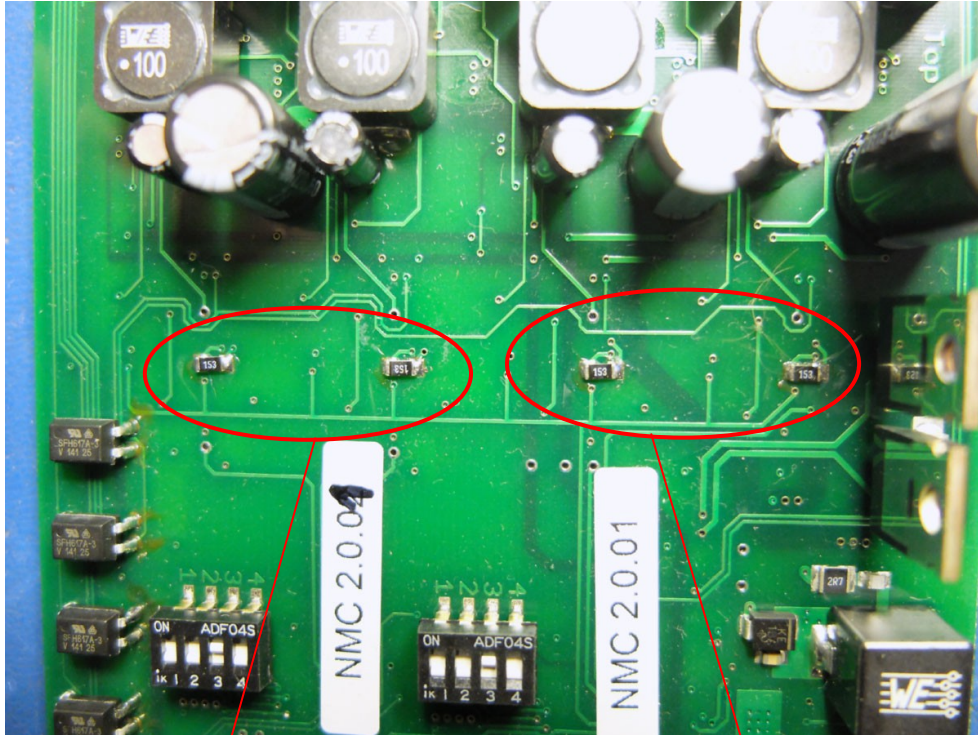

Current Resistors
(2) Channel 2

Current Resistors
(2) Channel 1

Resistor	Current
33k	0.96A
22k	1.46A
15k	2.15A
10k	3.24A
8.2k	3.87A

Resistors must be matched / paired within each axis!