

<b>Title:</b>	<b>DDC Error Codes</b>			<b>ID:</b>
				140
<b>22.10.14</b>	<b>Response:</b>	<b>Model:</b>	<b>Author:</b>	
		-	CMA	

## DDC Error Codes

Code	Error	Comment
<b>E01</b>	No Sensor	No response of sensor over a long period (10..20s) of time
<b>E02</b>	RX Timeout	No response on data-requests to sensor (1s)
<b>E03</b>	TX Timeout	No start-bit within time out
<b>E04</b>	NaN Error	Number parsing error
<b>E05</b>	Checksum Error	Check sum error over data string package

CMA, 22.10.14