

|                 |  |               |                |
|-----------------|--|---------------|----------------|
| <b>Title:</b>   | <b>How to Verify Defrost Operation</b> | <b>ID:</b>    |                |
|                 |  | 0085          |                |
| <b>Date in:</b> | <b>Response:</b>                       | <b>Model:</b> | <b>Author:</b> |
| 2009-11-11      | 2009-11-11                             | -             | CMa            |

## How to Verify Defrost Operation

**Q:**

How do I check defrost cycle operation

**A:**

To find out:

- 1 blow air through drain tube. Watch if tubing blocked.
- 2 Defrost system, Watch for water coming through drain tube.
  - 2.1 If water is flowing through tube the flow should spot at the time when the defrost cycle stops
  - 2.2 If there is no water check if drain tube is free by blowing air through drain tube
- 3 Empty drain pan. Then turn system off 12..24 hours.
- 4 Monitor amount of water in drain pan.