

Title:	REMP SSS Temperature Alarm Definitions			ID:
				0091
Date in:	Response:	Model:	Author:	
2010-05-19	2010-05-19	-	C <i>Ma</i>	

## REMP SSS Temperature Alarm Definitions

- Alarm Output for Temperature out of band is 600
- Alarm Set point DM900, value:  $(t+500)*32$  [1/10°C]
- Band for temperature is DM130 [1/10 °C]
- Delay for Temp is T130 [1h]
- Alarm turns on after T130 time [1/10s]. As soon as temp. is within band.
- Reset Alarm: "RS T130" resets alarm for T130 (1 hour)